

## **IMMUNOFLUORESCENCE OF NEURONS**

- fix neurons with 4%PFA and 4% sucrose in PBS, 20 minutes at RT.
- Wash 3 x 2min with PBS
- Permeabilize with Triton X-100 0.1% for 5 minutes.
- Was 2x 2minutes with PBS
- Block with 0.5% BSA (filtered) for 20 minutes.
- Incubate primary Ab for 60 minutes
- Wash 3 x 2min with PBS
- Incubate with Secondary (alexa-labeld) Ab for 60 minutes
- Wash 3 x 2min with PBS
- Image.